

Density S4 4.1.1.1002 Release Notes

Part Number: 30239-RN

Rev: A

OCT-27-2021-4.1.1.1002

1. ADDED: Manual Export Options:
 - Add User Defined File Name.
 - Add Save Header as First Row check box.
 - Add Create New File Option.
2. CHANGED: Auto Export Options:
 - Don't grey out Overwrite File, Append to File radio buttons.
3. Changed: "Create file" to "Create New File"
4. ADDED: File Naming Options:
 - Add one more option Attribute to File Naming Options.
 - Default
 - Attribute
 - User Defined
5. Changed: Show current User Defined options when the user selects "Attribute"
6. Added: "User Defined File Name" - Input Box for "User Defined" option.
7. BUG: Do proper validation and show appropriate message when the user selects different combination of Export option and File Naming Option.
8. BUG: Records with no measurement data must not stop the export. If a measurement has no data - ignore and move on. Do not create a header for this measurement record just go to the next record
9. ADDED: Export of Multiple results from Calibration Verification and Calibration Adjustment
 - Checkboxes was added for each result.
 - "Export" button was added.
 - User can select multiple results and can export selected results to file.
10. ADDED: Add Calibration and Verification Report Print at the end of Calibration and Verification.
 - Print button was added in final prompt. Preview will be opened when the user presses this button. User can print the calibration result from preview.
11. BUG:Calibrating Air Water on Density Meter : System recommended reject when we perform calibration, covers bottom menu.